

Development and Characterization of Microfabricated Disposable Platinum Electrodes for High Performance Ion Chromatography

J. Cheng, P. Jandik, and C. Pohl, Dionex Corporation, Sunnyvale, CA, USA


INTRODUCTION

Pt Working Electrodes: Background Information


- Detection of a broad range of electro active groups
 - Hydroxy, methoxy, or amine groups in aliphatic, aromatic, and carbohydrate compounds (such as alcohols, antibiotics, phenolic antioxidants, catechols, and carbohydrates)
 - Amino thiols and other S-containing species
 - Formic acid and –OH containing organic acids
 - Inorganic Anions (such as SO_3^{2-} , S^{2-} , N_3^- and CN^-).
- Knitted PTFE reaction coil is used between the chromatographic column and the detection cell.
- Current methods with conventional Pt electrodes are not as popular as methods with Au electrodes
 - Poor reproducibility of published methods
 - Relatively difficult to develop a suitable waveform for achieving a stable detection response over a long period of time.

Why Disposable Pt Working Electrodes?

- Predictable stability of response
- Easy installation
 - Fast equilibration (minimal delay after installation)
 - Excellent electrode-to-electrode reproducibility
- No conditioning required
 - No hand polishing
 - No activation waveforms

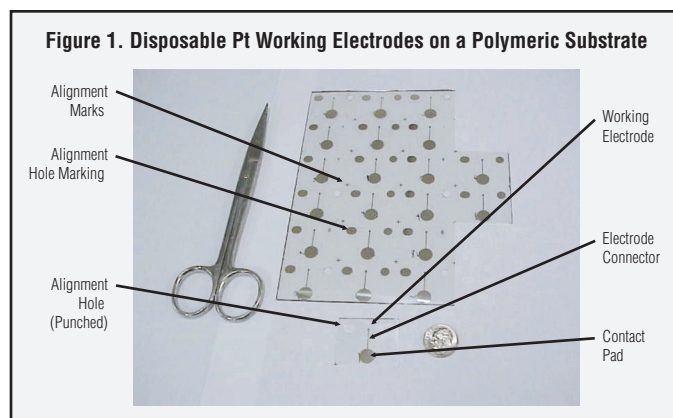

PULSED AMPEROMETRY WAVEFORM

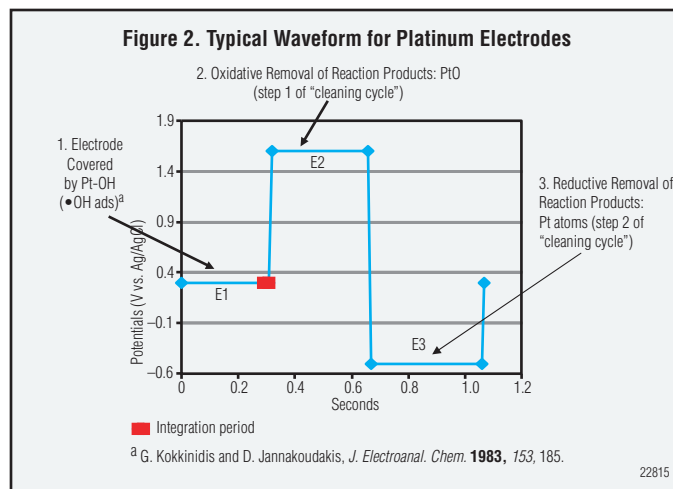


Now sold under the
Thermo Scientific brand

Thermo
SCIENTIFIC

NEW WAVEFORM DEVELOPMENT FOR DISPOSABLE PT ELECTRODES

Cyclic Voltammetry with Disposable Platinum Electrodes


Figure 3. Cyclic Voltammetry of Glycerol Performed Inside a Chromatographic Detection Cell of a Dionex ICS-3000

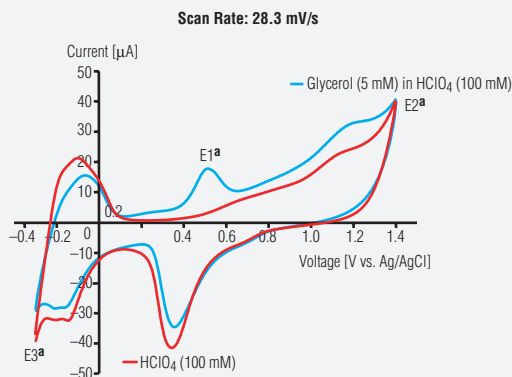

22816

Waveform Optimization

Figure 4. Detection Response vs. Detection Potential (E1)

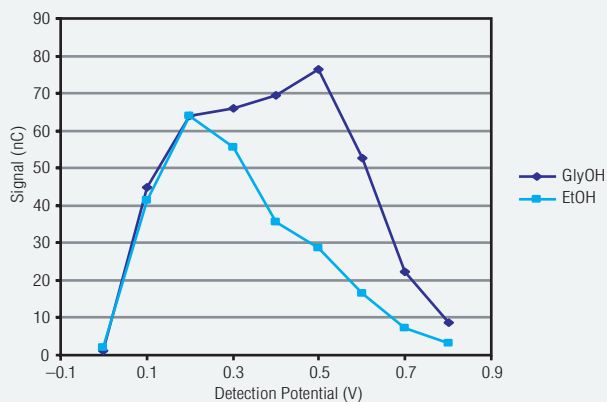

22817

Table 1. Effect of Oxidative Potential (E2 or Step 1 "Cleaning Cycle")

E2 ^a (V)	Peak Height (nC)		Peak Asymmetry		Lifetime ^b (Days)
	Glycerol	EtOH	Glycerol	EtOH	
1.00	43.10	12.15	1.83	2.52	11
1.10 ^c	45.40	12.91	1.86	1.96	11
1.20	47.02	13.92	2.09	1.76	-7
1.30	49.26	14.74	2.30	1.93	5-6
1.40	49.03	14.82	1.98	1.73	-3
1.50	50.05	16.56	1.83	1.17	N. D.
1.60	49.39	15.48	1.96	1.66	<1

^a In all runs, E1s and E3s are kept at 0.45 and -0.70 V, respectively.

^b The lifetime was defined as a time interval required to dissolve the platinum layer under chromatographic detection conditions.

^c Optimal conditions.

Table 2. Effect of Reductive Potential (E3 or Step 2 "Cleaning Cycle")


E3 ^a (V)	Peak Height (nC)		Peak Asymmetry		Noise (nC)
	Glycerol	EtOH	Glycerol	EtOH	
0.20	57.33	16.07	2.82	1.79	0.210
0.00	51.24	18.62	2.70	1.86	0.180
-0.20 ^b	47.71	18.54	2.48	1.52	0.147
-0.40 ^b	54.04	22.06	2.41	2.07	0.148
-0.50	52.37	22.66	2.42	2.07	0.151
-0.60	44.73	21.23	2.38	2.12	0.156
-0.80	44.55	22.73	2.51	2.11	0.143

^a Waveform conditions: detection potential (E1) is 0.45 V; Oxidation cleaning potential (E2) is 1.20 V.

^b Optimal conditions.

Figure 5. Varying Duration of Step 1 of the "Cleaning Cycle"

Concentrations: 50 ppm all analytes (except butanol: 100 ppm)

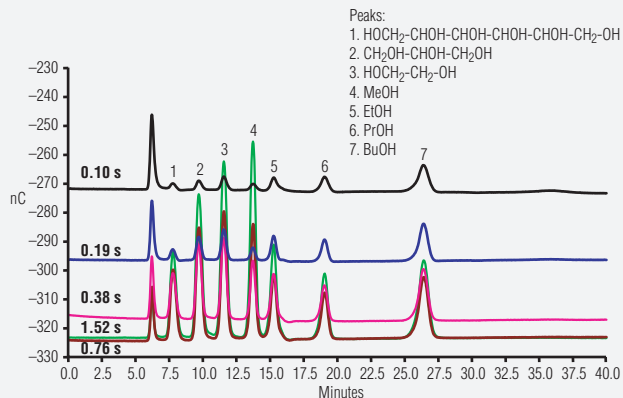

Cleaning Time (s)	Response (nC, GlyOH)	Ratio of S/N (GlyOH)
0.10	3	52
0.19	8	93
0.38	26	233
0.76	39	351
1.52	49	549

Total Cleaning Times: 1.52, 0.76, 0.38, 0.19 and 0.10 s

Total Waveform Times: 1.83, 1.07, 0.69, 0.50 and 0.41 s

22818

SEPARATION AND DETECTION METHODS




Time (s)	Potential (V) ^b	Integration
0.00	0.30	
0.29	0.30	Start
0.31	0.30	End
0.32	1.15	
0.66	1.15	
0.67	-0.30	
1.06	-0.30	
1.07	0.30	

^a RSD (% , electrode-to-electrode): Glycerol (3.46) and EtOH (12.35)

with the optimized waveform

^b Versus Ag/AgCl


ANALYTICAL PERFORMANCE OF DISPOSABLE PT ELECTRODES


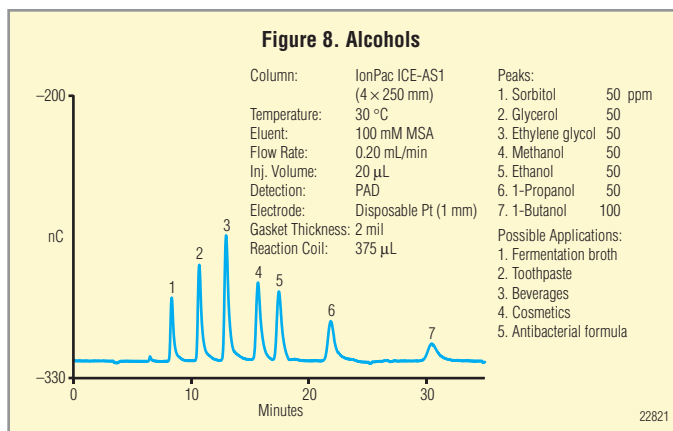
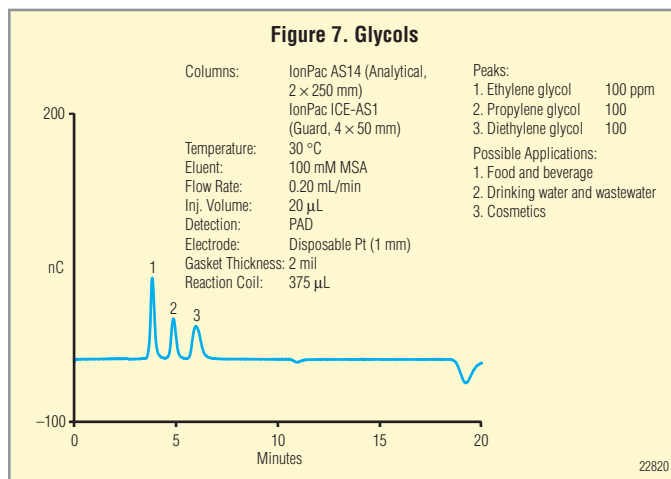
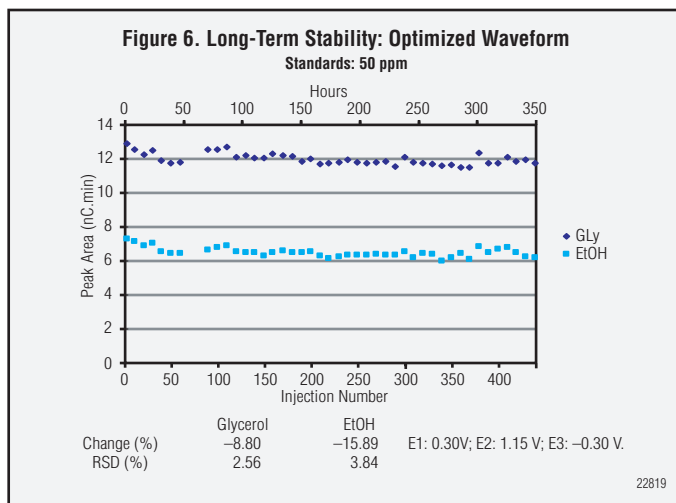


Alcohol	Electrode Type	Range (ppm)	Correlation Coefficient ^a
Glycerol	Non-disposable	0.5–100	1.0000
	Disposable	0.5–100	0.9995
Ethanol	Non-disposable	1–500	0.9996
	Disposable	1–500	0.9998


^a Quadratic fit

	Conventional Waveform	Optimized Waveform	
	ED 40 Pt (ppm)	ED 40 Pt (ppm)	Disposable Pt (ppm)
Alcohol			
Glycerol	0.54	0.20	0.10
Ethanol	1.63	0.51	0.25

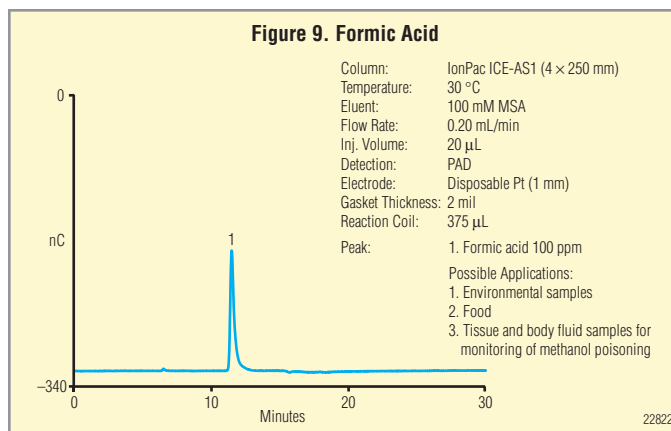

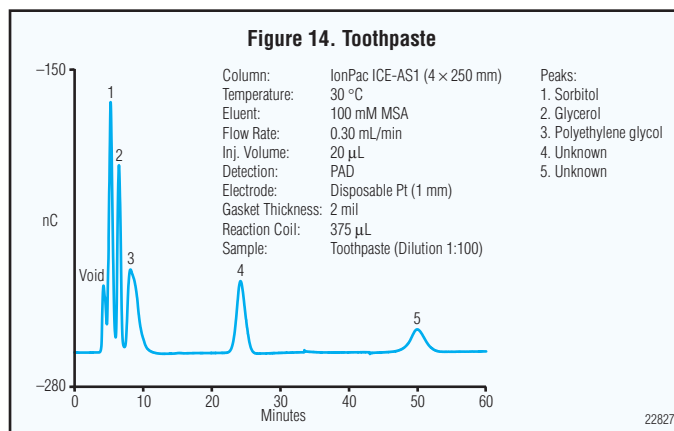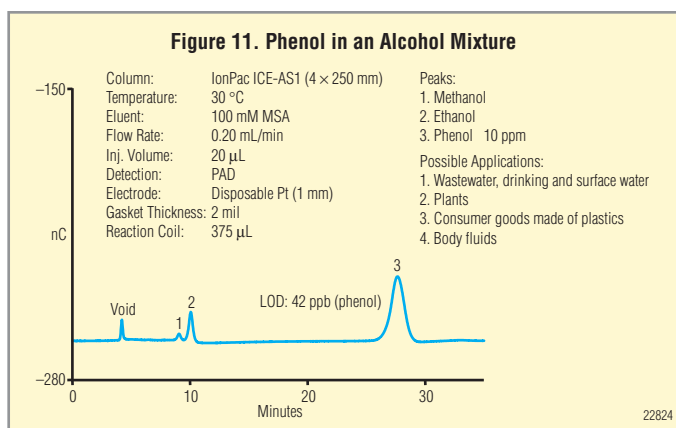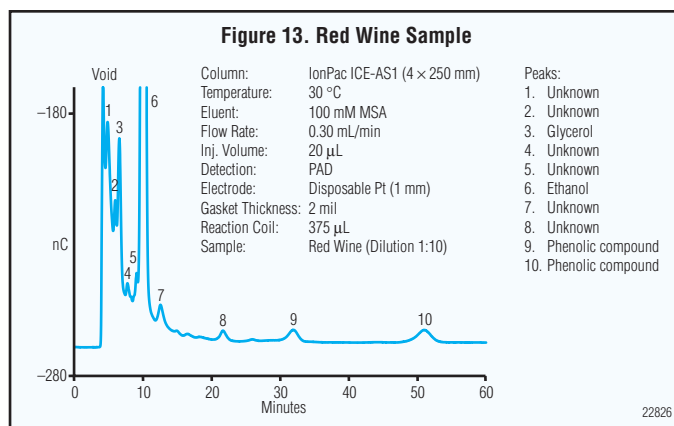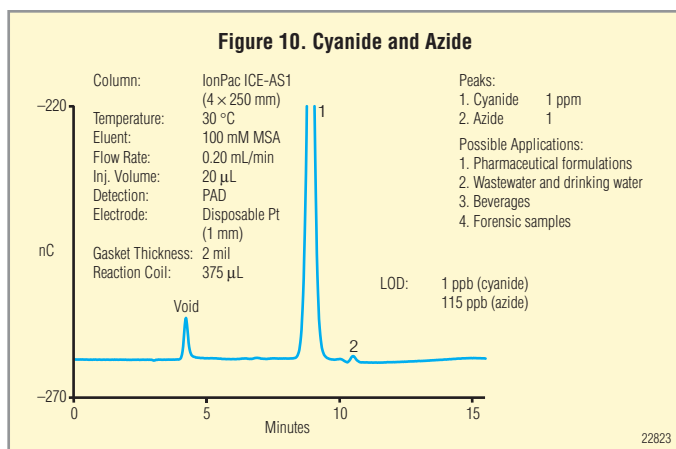

APPLICATIONS

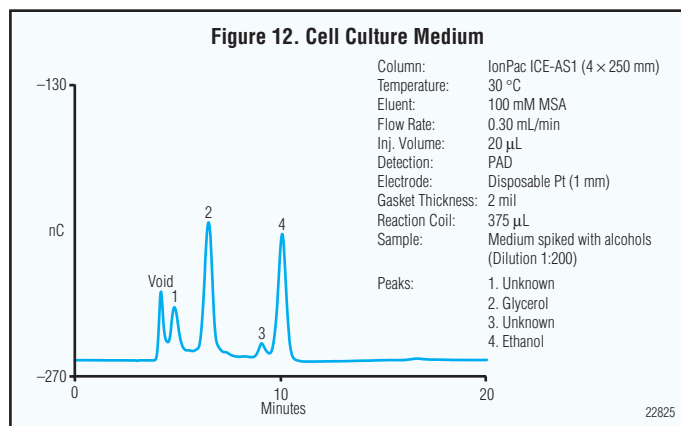

Conclusions

- Developed a fabrication process for manufacturing disposable Pt electrodes
 - Excellent electrode-to-electrode reproducibility
 - Faster start-up
 - No polishing and reconditioning
- Optimized a waveform for disposable and non-disposable Pt electrodes
- Improved analytical performance for determining alcohols and other compounds (e.g., cyanide and azide) in acidic eluent

IonPac is a registered trademark of Dionex Corporation.

Passion. Power. Productivity.



Dionex Corporation

1228 Titan Way
 P.O. Box 3603
 Sunnyvale, CA
 94088-3603
 (408) 737-0700

North America

Sunnyvale, CA (408) 737-8522 Westmont, IL (630) 789-3660
 Houston, TX (281) 847-5652 Atlanta, GA (770) 432-8100
 Marlton, NJ (856) 596-0600 Canada (905) 844-9650

www.dionex.com

Europe

Austria (43) 1 616 51 25 Belgium (32) 3 353 4294
 Denmark (45) 36 36 90 90 France (33) 1 39 30 01 10
 Germany (49) 6126 991 210 Italy (39) 06 66 51 5052
 The Netherlands (31) 161 43 43 03
 Switzerland (41) 62 205 99 66
 United Kingdom (44) 1276 691272

Asia Pacific

Australia (61) 2 9420 5233 China (852) 2428 3282
 India (91) 22 28475235 Japan (81) 6 6885 1213
 Korea (82) 2 2653 2580



LPN 1791-01 04/06
 ©2006 Dionex Corporation