

Easy Method Transfer from HPLC to RSLC with the Dionex Method Speed-Up Calculator

INTRODUCTION

The goal of every chromatographic optimization is a method that sufficiently resolves all peaks of interest in as short a time as possible. The evolution of packing materials and instrument performance has extended chromatographic separations to new limits: ultrahigh-performance liquid chromatography (UHPLC).

The new Dionex UltiMate® 3000 Rapid Separation LC (RSLC) system is ideal for ultrafast, high-resolution LC. The RSLC system was designed for ultrafast separations with flow rates up to 5 mL/min at pressures up to 800 bar (11,600 psi) for the entire flow-rate range. This industry-leading flow-pressure footprint ensures the highest flexibility possible; from conventional to ultrahigh-resolution to ultrahigh-speed methods. The RSLC system, with autosampler cycle times of only 15 seconds, oven temperatures up to 110 °C, and data

collection rates up to 100 Hz (even when acquiring UV-Vis spectra), sets the standard for UHPLC performance. Acclaim® RSLC columns with a 2.2 µm particle size complete the RSLC dimension.

A successful transfer from an HPLC method to an RSLC method requires recalculation of the chromatographic parameters. Underlying chromatographic principles have to be considered to find the appropriate parameters for a method transfer. With the Method Speed-up Calculator, Dionex offers an electronic tool that streamlines the process of optimum method transfer. This technical note describes the theory behind the Method Speed-Up Calculator and the application of this interactive, multi-language tool, illustrated with an exemplary method transfer from a conventional LC separation to an RSLC separation. You may obtain a copy of this calculator from your Dionex representative.

Now sold under the
Thermo Scientific brand

Thermo
SCIENTIFIC

METHOD SPEED-UP STRATEGY

The purpose of method speed-up is to achieve sufficient resolution in the shortest possible time. The strategy is to maintain the resolving power of the application by using shorter columns packed with smaller particles. The theory for this approach is based on chromatographic mechanisms, found in almost every chromatography text book. The following fundamental chromatographic equations are applied by the Method Speed-Up Calculator for the method transfer from conventional to ultrafast methods.

The separation efficiency of a method is stated by the peak capacity P , which describes the number of peaks that can be resolved in a given time period. The peak capacity is defined by the run time divided by the average peak width. Hence, a small peak width is essential for a fast method with high separation efficiency. The peak width is proportional to the inverse square root of the number of theoretical plates N generated by the column. Taking into account the length of the column, its efficiency can also be expressed by the height equivalent to a theoretical plate H . The relationship between plate height H and plate number N of a column with the length L is given by Formula 1.

$$\text{Formula 1: } N = \frac{L}{H}$$

Low height equivalents will therefore generate a high number of theoretical plates, and hence small peak width for high peak capacity is gained. Which factors define H ? For an answer, the processes inside the column have to be considered, which are expressed by the Van Deemter equation (Formula 2).

$$\text{Formula 2: } H = A + \frac{B}{u} + C \cdot u$$

The Eddy diffusion A describes the mobile phase movement along different random paths through the stationary phase, resulting in broadening of the analyte band. The longitudinal diffusion of the analyte against the flow rate is expressed by the term B . Term C describes the resistance of the analyte to mass transfer into the pores of the stationary phase. This results in higher band broadening with increasing velocity of the mobile phase. The well-known Van Deemter plots of plate height H against the linear velocity of the mobile phase are useful

in determining the optimum mobile phase flow rate for highest column efficiency with lowest plate heights. A simplification of the Van Deemter equation, according to Halász¹ (Formula 3), describes the relationship between column efficiency (expressed in plate height H), particle size d_p (in μm) and velocity of mobile phase u (in mm/s):

$$\text{Formula 3: } H = 2 \cdot d_p + \frac{6}{u} + \frac{d_p^2 \cdot u}{20}$$

The plots of plate height H against velocity u depending on the particle sizes d_p of the stationary phase (see Figure 1, top) demonstrate visually the key function of small particle sizes in the method speed-up strategy: The smaller the particles, the smaller the plate height and therefore the better the separation efficiency. An efficiency equivalent to larger particle columns can be achieved by using shorter columns and therefore shorter run times.

Another benefit with use of smaller particles is shown for the 2 μm particles in Figure 1: Due to improved mass transfer with small particle packings, further acceleration of mobile phases beyond the optimal flow rate with minimal change in the plate height is possible.

Optimum flow rates and minimum achievable plate heights can be calculated by setting the first derivative of the Halász equation to zero. The optimal linear velocity (in mm/s) is then calculated by Formula 4.

$$\text{Formula 4: } u_{opt} = \sqrt{\frac{B}{C}} = \frac{10.95}{d_p}$$

The minimum achievable plate height as a function of particle size is calculated by insertion of Formula 4 in Formula 3, resulting in Formula 5.

$$\text{Formula 5: } H_{min} \approx 3 \cdot d_p$$

Chromatographers typically prefer resolution over theoretical plates as a measure of the separation quality. The achievable resolution R of a method is directly proportional to the square root of the theoretical plate number as can be seen in Formula 6. k is the retention factor of the analyte and α the selectivity.

$$\text{Formula 6: } R = \frac{1}{4} \cdot \sqrt{N} \cdot \frac{k_2}{1+k_2} \cdot \frac{\alpha-1}{\alpha}$$

If the column length is kept constant and the particle size is decreased, the resolution of the analytes improves. Figure 1, bottom, demonstrates this effect using 5 μm and 2 μm particles.

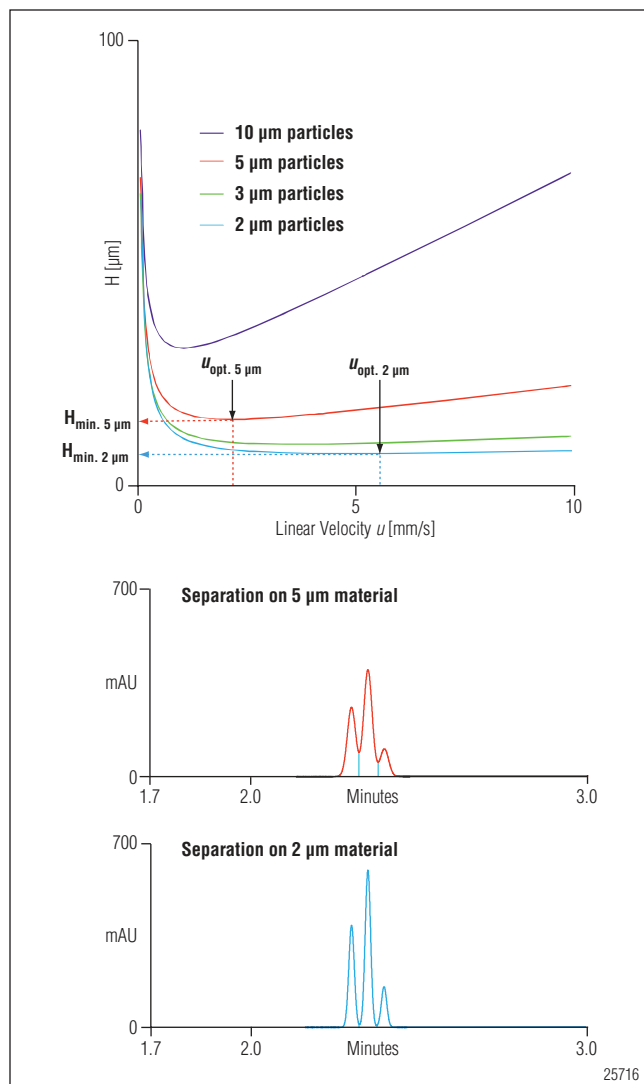


Figure 1. Smaller particles provide more theoretical plates and more resolution, demonstrated by the improved separation of three peaks (bottom) and smaller minimum plate heights H in the Van Deemter plot (top). At linear velocities higher than u_{opt} , H increases more slowly when using smaller particles, allowing higher flow rates and therefore faster separations while keeping separation efficiency almost constant. The speed-up potential of small particles is revealed by the Van Deemter plots (top) of plate height H against linear velocity u of mobile phase: Reducing the particle size allows higher flow rates and shorter columns because of the decreased minimum plate height and increased optimum velocity. Consequently, smaller peak width and improved resolution are the result (bottom).

When transferring a gradient method, the scaling of the gradient profile to the new column format and flow rate has to be considered to maintain the separation performance. The theoretical background was introduced by L. Snyder² and is known as the gradient volume principle. The gradient volume is defined as the mobile phase volume that flows through the column at a defined gradient time t_G . Analytes are considered to elute at constant eluent composition. Keeping the ratio between the gradient volume and the column volume constant therefore results in a correct gradient transfer to a different column format.

Taking into account the changed flow rates F and column volume (with diameter d_c and length L), the gradient time intervals t_G of the new methods are calculated with Formula 7.

$$\text{Formula 7: } t_{G,\text{new}} = t_{G,\text{old}} \cdot \frac{F_{\text{old}}}{F_{\text{new}}} \cdot \frac{L_{\text{new}}}{L_{\text{old}}} \cdot \left(\frac{d_{c,\text{new}}}{d_{c,\text{old}}} \right)^2$$

An easy transfer of method parameters can be achieved by using the Dionex Method Speed-Up Calculator (Figure 2), which incorporates all the overwhelming theory and makes manual calculations unnecessary. This technical note describes the easy method transfer of an example separation applying the calculator. Just some prerequisites described in the following section have to be taken into account.

PREREQUISITES

The Method Speed-Up Calculator is a universal tool and not specific for Dionex products. Nevertheless, some prerequisites have to be considered for a successful method transfer, which is demonstrated in this technical note by the separation of seven soft drink additives.

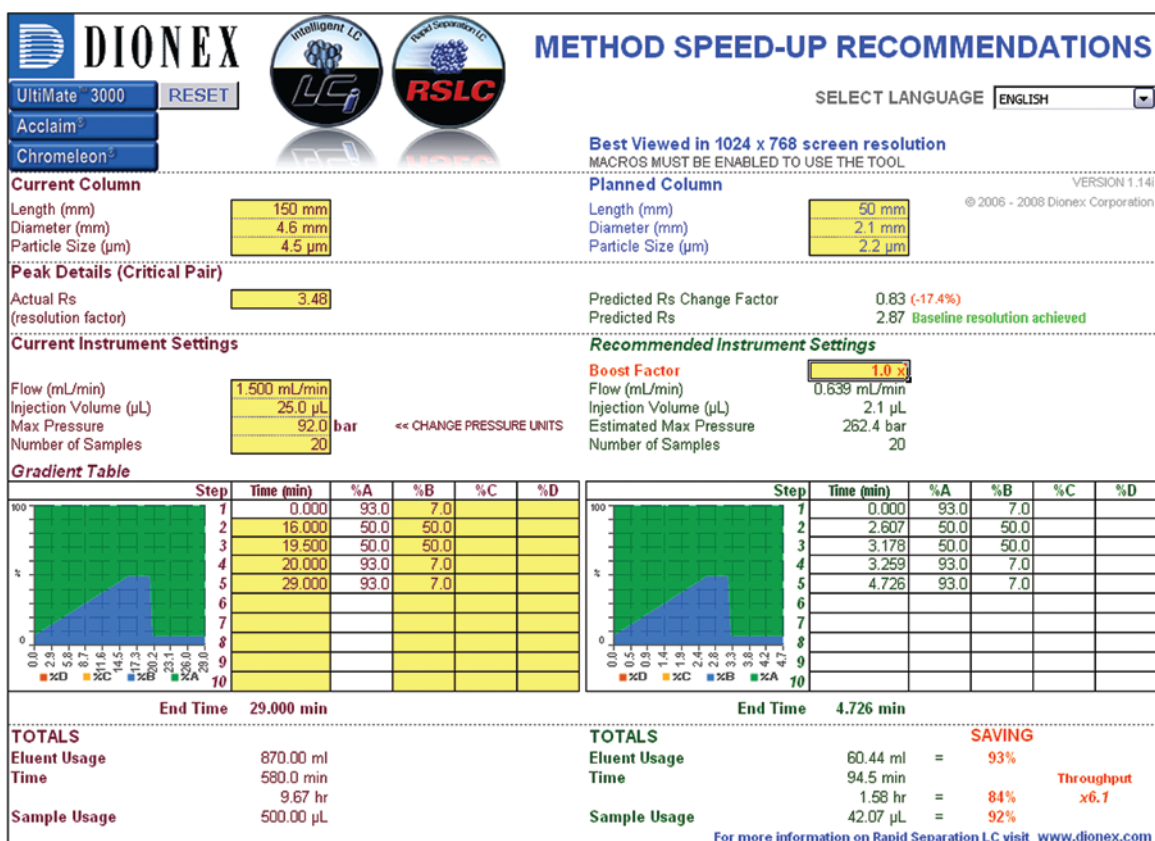


Figure 2. The Dionex Method Speed-Up Calculator transfers a conventional (current) HPLC method to a new (planned) RSLC method.

Column Dimension

First, the transfer of a conventional method to an RSLC method requires the selection of an adequate column filled with smaller particles. The RSLC method is predicted best if the selectivity of the stationary phase is maintained. Therefore, a column from the same manufacturer and with nominally identical surface modification is favoured for an exact method transfer. If this is not possible, a column with the same nominal stationary phase is the best choice. The separation is made faster by using shorter columns, but the column should still offer sufficient column efficiency to allow at least a baseline separation of analytes. Table 1 gives an overview of the theoretical plates expected by different column length and particle diameter size combinations using Dionex Acclaim column particle sizes. Note that column manufacturers typically fill columns designated 5 μm with particle sizes 4-5 μm. Dionex Acclaim 5 μm columns are actually filled with 4.5 μm particles. This is reflected in the table.

Table 1. Theoretical Plates Depending on Column Length and Particle Diameter (Calculated Using Formula 5)

	Theoretical Plates N		
	4.5 μm	3 μm	2.2 μm
Particle size	4.5 μm	3 μm	2.2 μm
Column length: 250 mm	18518	27778	37879
150 mm	11111	16667	22727
100 mm	7407	11111	15152
75 mm	5555	8333	11364
50 mm	3703	5556	7576

If the resolution of the original separation is higher than required, columns can be shortened. Keeping the column length constant while using smaller particles improves the resolution. Reducing the column diameter does not shorten the analysis time but decreases mobile phase consumption and sample volume. Taking into account an elevated temperature, smaller column inner diameters reduce the risk of thermal mismatch.

System Requirements

Smaller particles generate higher backpressure. The linear velocity of the mobile phase has to be increased while decreasing the particle size to work within the Van Deemter optimum. The UltiMate 3000 RSLC system perfectly supports this approach with its high maximum operation pressure of 800 bar (11,600 psi). This maximum pressure is constant over the entire flow rate range of up to 5 mL/min, providing additional potential to speed up applications even further by increasing the flow rate.

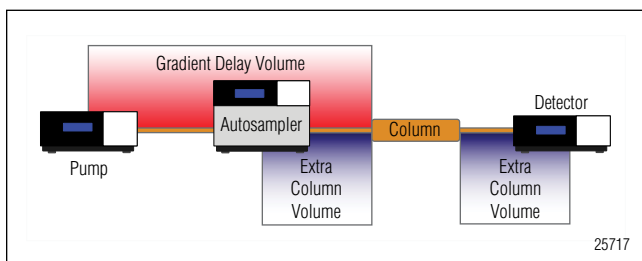


Figure 3. Gradient delay volume and extra column volume of an HPLC system. Both play an important role in method speed-up.

For fast gradient methods, the gradient delay volume (GDV) plays a crucial role. The GDV is defined as the volume between the first point of mixing and the head of the column. The GDV becomes increasingly important with fast, steep gradients and low flow rate applications as it affects the time taken for the gradient to reach the head of the column. The larger the GDV, the longer the initial isocratic hold at the beginning of the separation. Typically, this leads to later peak elution times than calculated. Early eluting peaks are affected most. In addition, the GDV increases the time needed for the equilibration time at the end of a sample and therefore increases the total cycle time. A general rule is to keep the gradient steepness and the ratio of GDV to column volume constant when transferring a standard method into a fast LC method. This will maintain the selectivity of the original method.³

The GDV can be adjusted to the column volume by installing appropriate mixer kits to the RSLC pump (see Table 2), which contributes most to the GDV. Typically, 100 μL or 200 μL mixers are good starting points when operating a small volume column in an RSLC system.

Table 2. Mixer Kits Available for UltiMate 3000 RSLC System to Adapt GDV of Pump

Mixer Kit	GDV pump
Mixer kit 6040.5000	35 μL
Static mixer kit 6040.5100	100 μL
Static mixer kit 6040.5150	200 μL

Another option is to switch the sample loop of the split-loop autosampler out of the flow path. The GDV is then reduced by the sample loop volume in the so-called bypass mode. The GDV of a standard sample loop of the RSLC autosampler is 150 μL , the micro injection loop has a 50 μL GDV.

Besides the gradient delay volume, the extra column volume is an important parameter for fast LC methods. The extra column volume is the volume in the system through which the sample passes and hence contributes to the band broadening of the analyte peak (Figure 3). The extra column volume of an optimized LC system should be below $1/10^{\text{th}}$ of the peak volume. Therefore the length and inner diameter of the tubing connections from injector to column and column to detector should be as small as possible. Special care has to be taken while installing the fittings to avoid dead volumes. In addition, the volume of the flow cell has to be adapted to the peak volumes eluting from the RSLC column. If possible, the flow cell detection volume should not exceed $1/10^{\text{th}}$ of the peak volume.

Detector Settings

When transferring a conventional method to an RSLC method, the detector settings have a significant impact on the detector performance. The data collection rate and time constant have to be adapted to the narrower peak shapes. In general, each peak should be defined by at least 30 data points. The data collection rate and time constant settings are typically interrelated to optimize the amount of data points per peak and reduce short-term noise while still maintaining peak height, symmetry, and resolution.

The Chromeleon[®] Chromatography Management Software has a wizard to automatically calculate the best settings, based on the input of the minimum peak width at half height of the chromatogram. This width is best determined by running the application once at maximum data rate and shortest time constant. The obtained peak width may then be entered into the wizard for optimization of the detection settings. Refer to the detector operation manual for further details.

METHOD SPEED-UP USING THE CALCULATOR

Separation Example

Separation was performed on an UltiMate 3000 RSLC system consisting of a HPG-3200RS Binary Rapid Separation Pump, a WPS-3000RS Rapid Separation Well Plate Sampler with analytical sample loop (100 μ L), a TCC-3000RS Rapid Separation Thermostatted Column Compartment with precolumn heater (2 μ L), and a VWD-3400RS Variable Wavelength Detector with semi-micro flow cell (2.5 μ L). Chromeleon Chromatography Management Software (version 6.80, SR5) was used for both controlling the instrument and reporting the data. The modules were connected with stainless steel micro capillaries, 0.01" ID, 1/16" OD when applying the conventional LC method, 0.007" and 0.005" ID, 1/16" OD when applying the RSLC methods. A standard mixture of seven common soft drink additives was separated by gradient elution at 45 °C on two different columns:

- Conventional HPLC Column: Acclaim 120, C18, 5 μ m, 4.6 \times 150 mm column, (P/N 059148)
- Rapid Separation Column: Acclaim RSLC 120, C18, 2.2 μ m, 2.1 \times 50 mm column (P/N 068981).

The UV absorbance wavelength at 210 nm was recorded at 5 Hz using the 4.6 \times 150 mm column and at 25 Hz and 50 Hz using the 2.1 \times 50 mm column. Further method details such as flow rate, injection volume, and gradient table of conventional and RSLC methods are described in the following section. The parameters for the method transfer were calculated with the Dionex Method Speed-Up Calculator (version 1.14i).

The conventional separation of seven soft drink additives is shown in Figure 4A. With the Method Speed-Up Calculator, the method was transferred successfully to RSLC methods (Figure 4B and C) at two different flow rates. The easy method transfer with this universal tool is described below.

Column Selection for Appropriate Resolution

The column for method speed-up must provide sufficient efficiency to resolve the most critical pairs. In this example, separating peaks 5 and 6 is most challenging. A first selection of the planned column dimensions can be made by considering the theoretical plates according to Table 1. The 4.6 \times 150 mm, 5 μ m column is actually filled with 4.5 μ m particles. Therefore, it provides 11,111 theoretical plates. On this column, the

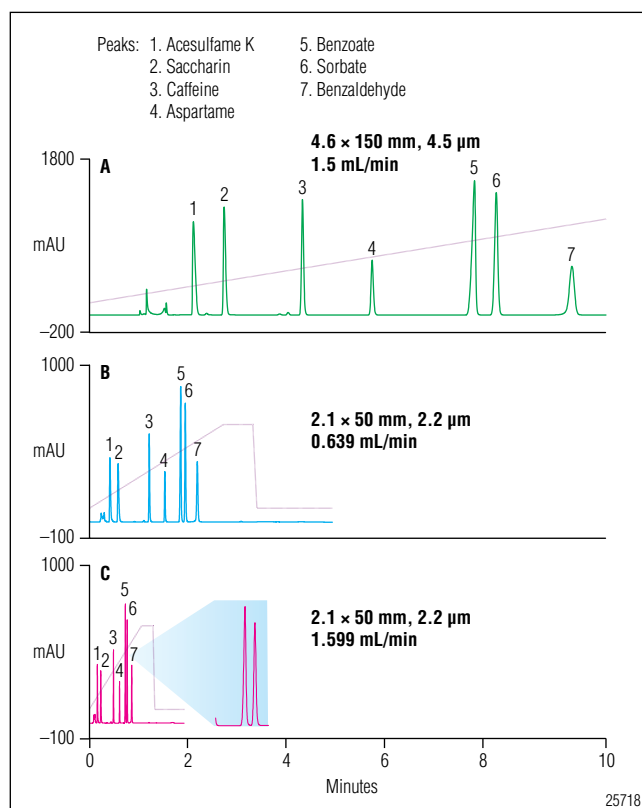


Figure 4. Method transfer with the Method Speed-Up Calculator from A) a conventional LC separation on an Acclaim 5 μ m particle column, to B) and C) RSLC separations on an Acclaim 2.2 μ m particle column.

resolution is $R(5,6)=3.48$. This resolution is sufficiently high to select a fast LC column with fewer theoretical plates for the speed up. Therefore, a 2.1 \times 50 mm, 2.2 μ m column with 7579 plates was selected.

The first values to be entered into the yellow field of the Method Speed-Up Calculator are the current column dimension, planned column dimension, and the resolution of the critical pair. To obtain the most accurate method transfer, use the particle sizes listed in the manufacturer's column specifications sheet instead of the nominal size, which may be different. Dionex Acclaim columns with a nominal particle size of 5 μ m are actually filled with 4.5 μ m particles, and this value should be used to achieve a precise method transfer calculation. This has a positive impact on the performance and pressure predictions for the planned column. Based on the assumption of unchanged stationary phase chemistry, the calculator then predicts the resolution provided by the new method (Figure 5).

Current Column		Planned Column		VERSION 1.14i © 2006 - 2008 Dionex Corporation
Length (mm)	150 mm	Length (mm)	50 mm	
Diameter (mm)	4.6 mm	Diameter (mm)	2.1 mm	
Particle Size (µm)	4.5 µm	Particle Size (µm)	2.2 µm	
Peak Details (Critical Pair)				
Actual Rs (resolution factor)	3.48	Predicted Rs Change Factor	0.83 (-17.4%)	
		Predicted Rs	2.87	Baseline resolution achieved

Figure 5. Column selection considering the resolution of the critical pair:

Current Instrument Settings		Recommended Instrument Settings	
Flow (mL/min)	1.500 mL/min	Boost Factor	1.0 x
Injection Volume (µL)	25.0 µL	Flow (mL/min)	0.639 mL/min
Max Pressure	92.0 bar	Injection Volume (µL)	2.1 µL
Number of Samples	20	Estimated Max Pressure	262.4 bar
	<< CHANGE PRESSURE UNITS	Number of Samples	20

Figure 6. The flow rate, injection volume and backpressure of the current method are scaled to the new column dimension.

In the example in Figure 5, the predicted resolution between benzoate and sorbate is 2.87. With a resolution of $R \geq 1.5$, the message “Baseline resolution achieved” pops up. This indicates that a successful method transfer with enough resolution is possible with the planned column. If R is smaller than 1.5, the red warning “Baseline is not resolved” appears. Note that the resolution calculation is performed only if the boost factor BF is 1, otherwise it is disabled. The function of the boost factor is described in the Adjust Flow Rate section.

Instrument Settings

The next section of the Method Speed-Up Calculator considers basic instrument settings. These are flow rate, injection volume, and system backpressure of the current method (Figure 6). In addition to these values, the detector settings have to be considered as described in the earlier section “Detector Settings”. Furthermore, the throughput gain with the new method can be calculated if the number of samples to be run is entered.

Adjust Flow Rate

As explained by Van Deemter theory, smaller particle phases need higher linear velocities to provide optimal separation efficiency. Consequently, the Dionex Method Speed-Up Calculator automatically optimizes the linear velocity by the ratio of particle sizes of the current and planned method. In addition, the new flow rate is scaled to the change of column cross section if the column

inner diameter changed. This keeps the linear velocity of the mobile phase constant. A boost factor (BF) can be entered to multiply the flow rate for a further decrease in separation time. If the calculated resolution with $BF=1$ predicts sufficient separation, the method can be accelerated by increasing the boost factor and therefore increasing the flow rate. Figure 1 shows that applying linear velocities beyond the optimum is no problem with smaller particle phases, as they do not significantly loose plates in this region. Note that the resolution calculation of the Method Speed-Up Calculator is disabled for $BF \neq 1$.

For the separation at hand, the flow rate is scaled from 1.5 mL/min to 0.639 mL/min when changing from an Acclaim 4.6 × 150 mm, 4.5 µm column to a 2.1 × 50 mm, 2.2 µm column (see Figure 6), adapting the linear velocity to the column dimensions and the particle size. The predicted resolution between peak 5 and 6 for the planned column is $R=2.87$. The actual resolution achieved is $R=2.91$, almost as calculated (chromatogram B in Figure 4).

A Boost Factor of 2.5 was entered for further acceleration of the method (Figure 7). The method was then performed with a flow rate of 1.599 mL/min, and resolution of the critical pair was still sufficient at $R=2.56$ (see zoom in chromatogram C in Figure 4).

Current Instrument Settings		Recommended Instrument Settings	
Flow (mL/min)	1.500 mL/min	Boost Factor	2.5 x 0.639 mL/min
Injection Volume (µL)	25.0 µL	Flow (mL/min)	1.599 mL/min
Max Pressure	92.0 bar	Injection Volume (µL)	2.1 µL
Number of Samples	20	Estimated Max Pressure	656.1 bar
Number of Samples	20	Number of Samples	20

<< CHANGE PRESSURE UNITS

Figure 7. The new flow rate is further accelerated by applying the Boost Factor of 2.5.

Scale Injection Volume

The injection volume has to be adapted to the new column dimension to achieve similar peak heights by equivalent mass loading. Therefore the injection plug has to be scaled to the change of column cross section. In addition, shorter columns with smaller particles cause a reduced zone dilution. Consequently, sharper peaks compared to longer columns are expected. The new injection volume $V_{inj,new}$ is then calculated by Formula 8, taking a changed cross section and reduced band broadening by changed particle diameter into account.

$$\text{Formula 8: } V_{inj,new} = V_{inj,old} \cdot \left(\frac{d_{c,new}}{d_{c,old}} \right)^2 \cdot \sqrt{\frac{L_{new} \cdot d_{p,new}}{L_{old} \cdot d_{p,old}}}$$

Generally, it is recommended that a smaller flow cell be used with the RSLC method to minimize the extra column volume. Also, the difference in path length of different flow cell sizes has to be taken into account while scaling the injection volume. In the example of the soft drink analysis, the injection volume is scaled from 25 µL to 2.1 µL when replacing the Acclaim 4.6 × 150 mm, 4.5 µm column with a 2.1 × 50 mm, 2.2 µm column (see Figure 6).

Predicted Backpressure

Speeding-up the current method by decreasing particle size and column diameter and increasing flow rate means elevating the maximum generated backpressure. The pressure drop across a column can be approximated by the Kozeny-Carman formula.⁴ The pressure drop of the new method is predicted by the calculator considering changes in column cross section, flow rate, and particle size and is multiplied by the boost factor. The viscosity of mobile phase is considered constant during method

transfer. The calculated pressure is only an approximation and does not take into account nominal and actual particle size distribution depending on column manufacturer. If the predicted maximum pressure is above 800 bar (11,600 psi) the warning “Exceeds pressure limit RSLC” is shown, indicating the upper pressure limit of the UltiMate 3000 RSLC system. However, in the case the method is transferred to a third party system, its pressure specification has to be considered.

In the example of the soft drink analysis, the actual pressure increases from 92 bar to 182 bar with $BF=1$ on the 2.1 × 50 mm column, and to 460 bar for the RSLC method with $BF=2.5$. The pressures predicted by the Method Speed-Up Calculator are 262 bar and 656 bar, respectively. The pressure calculation takes into account the change of the size of the column packing material. In a speed up situation, the pressure is also influenced by other factors such as particle size distribution, system fluidics pressure, change of flow cell, etc. When multiplication factors such as the boost factor are used, the difference between calculated and real pressure is pronounced. The pressure calculation is meant to give an orientation, what flow rates might be feasible on the planned column. However, it should be confirmed by applying the flow on the column.

Adapt Gradient Table

The gradient profile has to be adapted to the changed column dimensions and flow rate following the gradient-volume principle. The gradient steps of the current method are entered into the yellow fields of the gradient table. The calculator then scales the gradient step intervals appropriately and creates the gradient table of the new method.

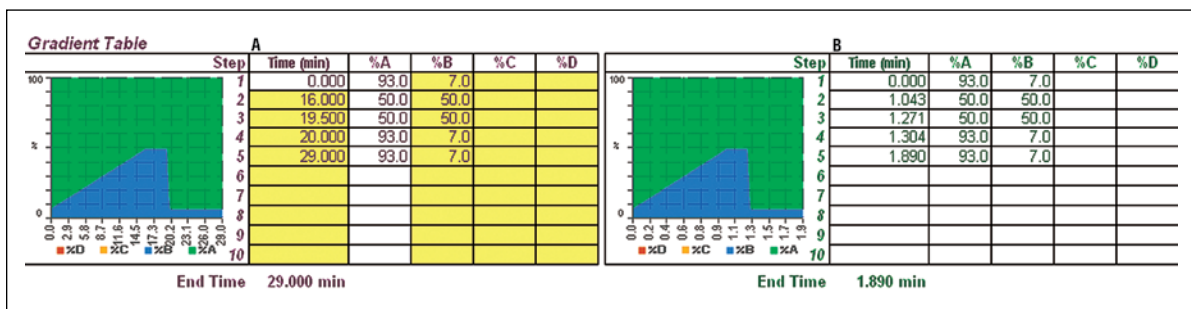


Figure 8. The gradient table of the current method (A) is adapted to the boosted method (B) according to the gradient-volume principle.

TOTALS					
Eluent Usage	870.00 ml	60.44 ml	=	93%	
Time	580.0 min	37.8 min	=	93%	Throughput
Sample Usage	500.00 μ L	42.07 μ L	=	92%	x15.3

Figure 9. The absolute values for analysis time, eluent usage, and sample usage of the current (purple) and planned (green) method are calculated by the Method Speed-Up Calculator. The savings of eluent, sample, and time due to the method transfer are highlighted.

The adapted gradient table for the soft drink analysis while using a boost factor $BF=1$ is shown in Figure 8. According to the gradient-volume principle, the total run time is reduced from 29.0 min to 4.95 min by taking into account the changed column volume from a 4.6×150 mm, $5 \mu\text{m}$ ($4.5 \mu\text{m}$ particles entered) to a 2.1×50 mm, $2.2 \mu\text{m}$ column and the flow rate reduction from 1.5 mL/min to 0.639 mL/min. The separation time was further reduced to 1.89 min by using boost factor $BF=2.5$. Gradient time steps were adapted accordingly. The comparison of the peak elution order displayed in Figure 4 shows that the separation performance of the gradient was maintained during method transfer.

Consumption and Savings

Why speed-up methods? To separate analyte peaks faster and at the same time reduce the mobile phase and sample volume consumption. Those three advantages of a method speed-up are indicated in the Method Speed-Up Calculator sheet right below the gradient table. The absolute values for the time, eluent, and sample usage are calculated taking the numbers of samples entered into the current instrument settings section of the calculation sheet into account (see Figure 6).

Regarding the soft drink analysis example, geometrical scaling of the method from the conventional column to the RSLC method means saving 93% of eluent and 92% of sample. The sample throughput increases 6.1-fold using $BF=1$. The higher flow rate at $BF=2.5$ results in a 15.3-fold increased throughput compared to the conventional LC method (Figure 9).

CONCLUSION

Fast method development or increased sample throughput are major challenges of most analytical laboratories. A systematic method speed-up is accomplished by reducing the particle size, shortening the column length, and increasing the linear velocity of the mobile phase. The Dionex Method Speed-Up Calculator automatically applies these rules and scales the conventional LC parameters to the conditions of the RSLC method. The interactive electronic tool is universally applicable. New instrument settings are predicted and gradient tables are adapted for optimum performance for the new method. The benefit of the method transfer is summarized by the integrated calculation of savings in time, eluent and sample. In addition, users can benefit from getting results earlier and thereby reducing the time to market. The Dionex Method Speed-Up Calculator is part of Dionex's total RSLC solution, which further consists of the industry leading UltiMate 3000 RSLC system, powerful Chromeleon Chromatography Management Software, and high-efficiency Acclaim RSLC columns.

REFERENCES

1. Halász, I.; Ende, R.; Asshauer, J. *J. Chromatogr., A* **1975**, *112*, 37–60.
2. Snyder, L.R.; Dolan, J.W.; Grant, J.R. *J. Chromatogr., A* **1979**, *165*, 3–30.
3. Schellinger, A.P.; Carr, P.W. *J. Chromatogr., A* **2005**, *1077*, 110-119.
4. Bear, J. *Dynamics of Fluids in Porous Media*; Dover: Mineola, NY, 1988.

Acclaim, Chromeleon, and UltiMate are registered trademarks of Dionex Corporation.

Passion. Power. Productivity.



Dionex Corporation

1228 Titan Way
P.O. Box 3603
Sunnyvale, CA
94088-3603
(408) 737-0700

North America

U.S. (847) 295-7500

South America

Brazil (55) 11 3731 5140

Europe

Austria (43) 1 616 51 25 Benelux (31) 20 683 9768 (32) 3 353 4294
Denmark (45) 36 36 90 90 France (33) 1 39 30 01 10 Germany (49) 6126 991 0
Ireland (353) 1 644 0064 Italy (39) 02 51 62 1267 Sweden (46) 8 473 3380
Switzerland (41) 62 205 9966 United Kingdom (44) 1276 691722

Asia Pacific

Australia (61) 2 9420 5233 China (852) 2428 3282 India (91) 22 2764 2735
Japan (81) 6 6885 1213 Korea (82) 2 2653 2580 Singapore (65) 6289 1190
Taiwan (886) 2 8751 6655

www.dionex.com



LPN 2140-01 PDF 2/09
©2009 Dionex Corporation