

Determination of Residual Acylglycerols in Biodiesel

INTRODUCTION

Many methods have been proposed to determine acylglycerols in biodiesel samples. This analysis is commonly performed by gas chromatography/flame ionization detection (GC/FID) methods such as ASTM D6584 and EN14105. However, GC methods often require sample derivatization, and high inlet temperatures can cause thermal degradation and underestimation of the sample amount. High-performance liquid chromatography (HPLC) analysis with a universal detector, such as charged aerosol detection (CAD), is emerging as the preferred method for biodiesel analysis. This technique is well suited to determine components with low volatility relative to the mobile phase used for separation.

Previous work has demonstrated that the ASTM GC method and HPLC methods can be statistically equivalent.¹ HPLC also permits direct analysis of biodiesel samples with just a simple dilution, precluding the derivatization step that is required for GC analysis of nonvolatile samples. Additionally, by using a small particle-size column, resolution can be improved and run times shortened.

This study demonstrates use of a rapid separation liquid chromatography (RSLC) method for analysis of acylglycerols in biodiesel fuel in under 10 min. This method avoids exceptionally toxic or environmentally damaging solvents, such as propionitrile and dichloromethane, which have been used for acylglycerol determination.²

Now sold under the
Thermo Scientific brand

Thermo
SCIENTIFIC

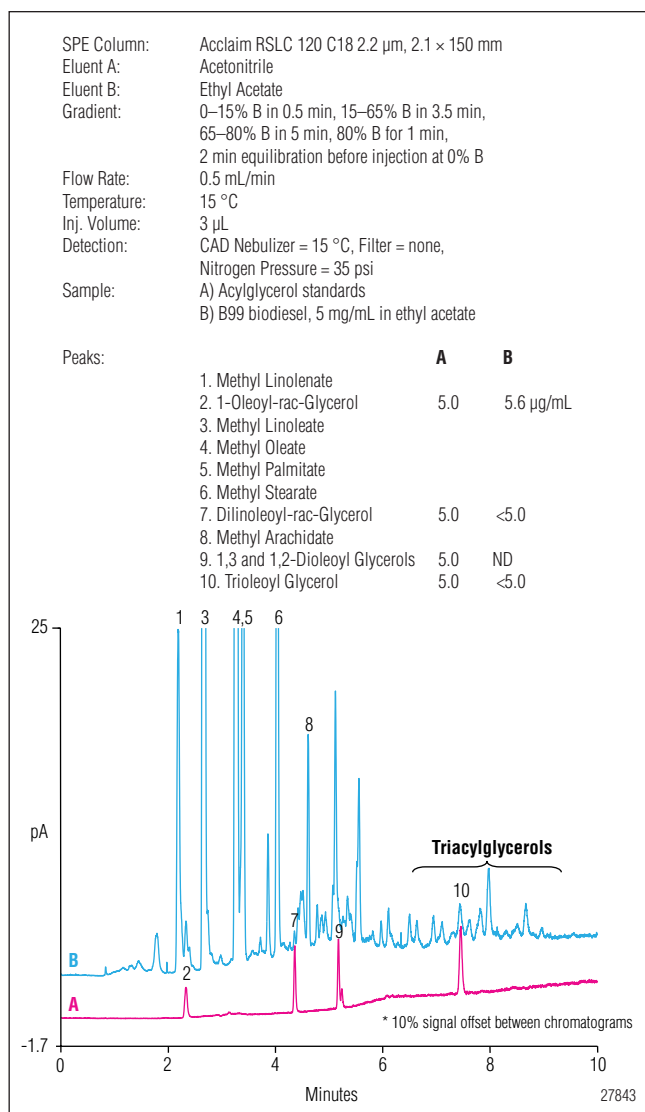


Figure 1. Separation of acylglycerols on the Acclaim RSLC 120 C18 column.

Figure 1 shows the separation of a B99 biodiesel sample produced from used fryer oil (B) and a mixture of acylglycerol standards (A). The triacylglycerols elute later than the fuel methyl esters and are well separated and easily detected using the Acclaim[®] RSLC 120 C18 column and the Corona[®] *ultra*[™] Charged Aerosol Detector (CAD). The wide dynamic range of the Corona CAD permits the detection of low-abundance oleoylglycerols in the

presence of highly concentrated components in a single analytical run. The use of UHPLC methods on the RSLC system with CAD detection has the sensitivity to determine acylglycerols in biodiesel samples at and below the 0.24% and 0.25% total glycerol limits in ASTM D6751 and EN 14214, respectively. This method provides a rapid and sensitive acylglycerols determination in biofuels with minimal sample preparation.

EQUIPMENT

System – UltiMate[®] RSLC with an HPG-3400RS pump, Corona *ultra* CAD, and WPS-3000TRS autosampler.

See Figure 1 for chromatography conditions.

Sample Preparation – Prepare biodiesel samples by dilution to 5 mg/mL in ethyl acetate prior to analysis.

REFERENCE

1. Foglia, T.A.; Jones, K.C.; Nunez, A.; Phillips, J.G.; and Mittelbach, M. Comparison of Chromatographic Methods for the Determination of Bound Glycerol in Biodiesel. *Chromatographia* **2004**, *60*, 305.
2. Buchgraber, M.; Ulberth, F.; Emons, H.; Anklam, E. Triacylglycerol Profiling by Using Chromatographic Techniques. *Eur. J. Lipid Sci. Technol.* **2004**, *106*, 621.

Acclaim, Corona, and UltiMate are registered trademarks and *ultra* is a trademark of Dionex.

Passion. Power. Productivity.



Dionex Corporation

1228 Titan Way
P.O. Box 3603
Sunnyvale, CA
94088-3603
(408) 737-0700

North America

U.S./Canada (847) 295-7500

South America

Brazil (55) 11 3731 5140

Europe

Austria (43) 1 616 51 25 Benelux (31) 20 683 9768 (32) 3 353 4294
Denmark (45) 36 36 90 90 France (33) 1 39 30 01 10 Germany (49) 6126 991 0
Ireland (353) 1 644 0064 Italy (39) 02 51 62 1267 Sweden (46) 8 473 3380
Switzerland (41) 62 205 9966 United Kingdom (44) 1276 691722

Asia Pacific

Australia (61) 2 9420 5233 China (852) 2428 3282 India (91) 22 2764 2735
Japan (81) 6 6885 1213 Korea (82) 2 2653 2580 Singapore (65) 6289 1190
Taiwan (886) 2 8751 6655

www.dionex.com



LPN 2626 PDF 10/10
©2010 Dionex Corporation